
1-800-295-5510 uline.com

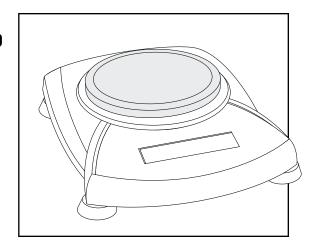

TROUBLESHOOTING

OPERATING ISSUE	RECOMMENDATIONS
No Display.	 Make sure the AC Adapter is plugged in. You may need to replace the batteries.
Battery Indicator is flashing.	Replace batteries.
Incorrect weight reading.	 Press the ON/ZERO OFF button with no weight on the platform, then weigh item. Recalibrate the balance.
Calibration procedures do not work.	 Use correct masses. See error codes note ERR4. (See Page 2)
Unable to display weight in a particular weighing unit.	 Use Units menu to set desired units ON. Exit Hold or Totalize mode.
Balance readings unstable.	Either move or shield the balance from external air currents or vibration.
Error code is displayed.	Review error code table and take appropriate action.

PAGE 1 OF 2 0113 TH-2382

uline.com

ERROR CODES

ERROR CODE LIST	RECOMMENDATIONS
Err 1 chEct	Cycle balance ON/OFF
Invalid checksum data.	2. Return balance for service
Err 2 LoAd	1. Remove load
Overload (>cap+90d) or Underload.	2. Make sure shipping lock is unlocked
	3. Recalibrate
Err 4 WElght	Use correct calibration weights
Incorrect calibration weight.	2. Check environmental conditions
Err 5 999999	Exit application
Displayed value>99999.	2. Rezero balance
Err 6 count	Exit application
Parts counting error – balance exits parts counting.	2. Rezero balance
Err 8 FULL	Cycle ON/OFF
RS232 buffer is full.	2. Check RS-232 settings in balance and computer
Err 9 dAtA	Cycle ON/OFF
Internal data error. The next button press or a 5 second time-out causes the balance to turn off.	2. Return balance for servicing